

INSTRUCTIONS FOR USE OF
CLASS II BIOLOGICAL SAFETY CABINET

TURN BLOWER ON 15 MINUTES PRIOR TO USE
DISINFECT INTERIOR SURFACES BEFORE WORK
COMMENCES

PLACE EQUIPMENT/SUPPLIES INTO CABINET

ALLOW 2-3 MINUTES FOR THE CABINET TO
PURGE AFTER PLACING MATERIALS IN IT

SLEEVES MUST BE WORN WHILE WORKING

DISINFECT INTERIOR SURFACES UPON
COMPLETION OF WORK

SC #044.2
Eff. 6/17

4 X 6