



6. To **start cycle**, press a numbered button at the right of the chamber. Press once to review preset times and temperatures. Press twice to start cycle.
7. The appropriate cycles for the corresponding loads are listed below:
  - a. Cycle 1: 65 min. @ 250<sup>0</sup>, dry- 25 min, soiled cages/ bedding.
  - b. Cycle 2: 30 min. @ 250<sup>0</sup>, dry- 15 min, clean cages & instruments.
  - c. Cycle 3: 30 min. @ 250<sup>0</sup>, dry- 25 min, double wrapped packs.
  - d. Cycle 4: 20 min. @ 250<sup>0</sup>, dry- 0 min, liquids, water bottles.
8. A beep will sound when the cycle is completed.
9. After the pressure in the chamber reaches zero the door can be opened by turning the handle counterclockwise and the load carefully removed. Vent the door by